## **Product Information**

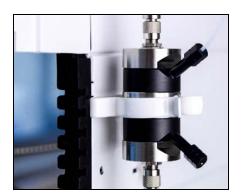


Fig. 1 PLATINblue flow cell mounted on detector side panel

The accessory kit for PLATINblue flow cells (Order no. F6911) with AZURA MWD 2.1L, DAD 2.1L and DAD 6.1L detectors enables the flow cell to be mounted onto the detector's housing via a spring clip and connected to the optics via fiber optic cables.

### Installation

#### Tools

Allen screwdriver size 2.5.

#### **Procedure**

- 1. Align the spring clip with the bore holes in the side panel.
- 2. Tighten the screw with the Allen screwdriver size 2.5.
- 3. Clip the flow cell into place.
- 4. Insert the fiber optics adapter.
- 5. Connect the fiber optic cables as shown in Fig.2.
- 6. Connect the capillaries.

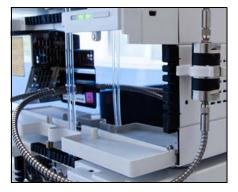


Fig. 2 Flow cell mounted on AZURA detector

# Part list of accessory kit F6911

Name	Order no.
Spring clip kit including clip and screw	A9855
Fiber optics cable 750 mm	A0742
Fiber optics cable 400 mm	A0911
Fiber optics adapter	AMKX8
Waste tubing kit	A9843
Capillaries with screw fittings	A2522 M2632 M2633

Note: For further details, consult the detector's user manual.