Read the corresponding technical documentation for handling and safety reasons.

Fitting guide for 50 mm ID columns

- Ensure to use the correct fittings for the 1/8" ports in the ID=50 mm columns.
- The ID=50 mm columns can only be connected with the corresponding UNF 5/16-24 screw threads.
- The usage of other fittings may lead to leakage and damage of the column inlets.
- Fittings for PEEK capillaries are included in the article A7226 (released for up to 100 bar).
- The fittings for stainless steel capillaries are included in the A7227 (released for up to 400 bar).



Any damage derived from the failure to follow these notes may void the columns warranty.

Fitting guide for 50 mm ID columns, V6759, version 2.0