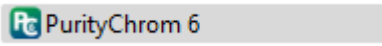
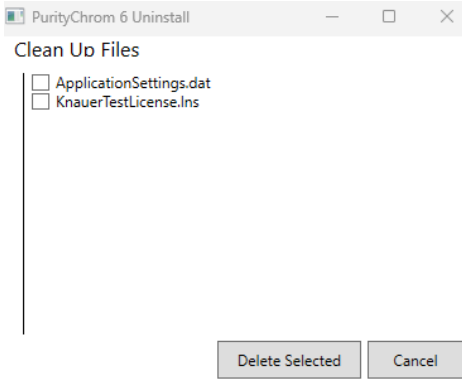
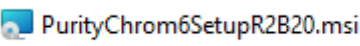
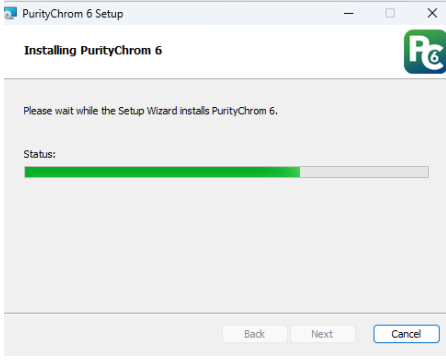


Updating the software version for PurityChrom 6

Note: All pictures of the software and software version numbers used in this instruction are exemplary and may vary in the PurityChrom 6 version you are using.

Process	Figure
<ol style="list-style-type: none"> 1. Open the Control Panel in Microsoft Windows. 2. Navigate to "Programs and Features". 	
<ol style="list-style-type: none"> 3. Locate and select „PurityChrom 6“ in the program list. 4. Right click on the selected icon and choose „Uninstall“. 5. The uninstallation wizard will start. 6. Confirm the dialog in the on-screen prompt. 	 <p>Fig. 1: Program icon of PurityChrom 6 in „Programs and features“</p>
<ol style="list-style-type: none"> 7. A subsequent dialog will appear, asking which files should be deleted. <p>Note: If you wish to maintain your license and configuration data, do not select any of the check boxes. Then click the „Delete Selected“ button. <ol style="list-style-type: none"> 8. The uninstallation process will be initiated. </p>	 <p>Fig. 2: Subsequent dialog</p>
<ol style="list-style-type: none"> 9. Install the latest version of PurityChrom 6 by either using the installer (provided on the software CD) or by executing the respective MSI file which is included in the delivery. 	 <p>Fig. 3: PurityChrom 6 installation file</p>
<ol style="list-style-type: none"> 10. Follow the on-screen prompts in the installation wizard in order to complete the process. 	 <p>Fig. 4: Installation dialog</p>

Process

Figure

11. Ensure that all installation steps (2 - 9) are marked „OK“.

Note: For a detailed description of each installation step, see the software manual ([V2670](#)).

12. Close the configuration wizard and click „Finish“ to complete the installation.

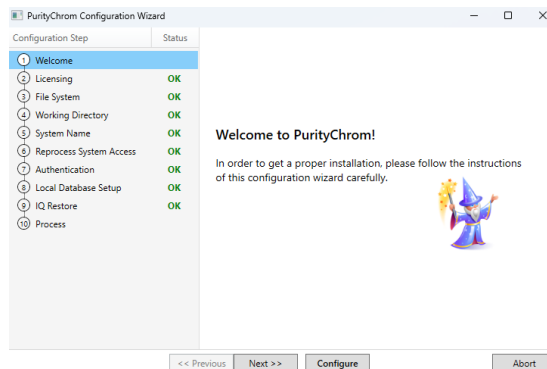



Fig. 5: Configuration wizard

13. Start PurityChrom 6.

14. To check if the installation has been successful, open the software settings  and check the displayed version number.

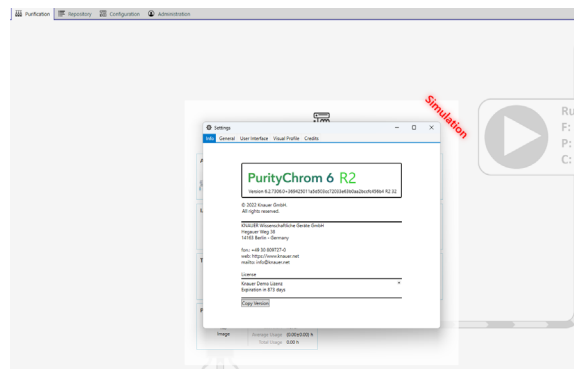


Fig. 6: Actual version number in the software

If you have followed the instruction, you will have uninstalled the previous PurityChrom 6 version and installed the new version successfully.